Contextual fear conditioning Procedure

Day 1. Conditioning training

Set up computer program and equipment (Acrylic chamber with shock grid clean it first using 70% of ethanol) for conditioning training. Set up light (100 Lux) using light meter. Set up shock 0.3miliampere and 65 dB for tone.

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1. Mouse handling for 1 minute then placed mouse carefully to the conditioning chamber.
2. Habituation of chamber before actual training for 120s.
3. Start the conditioning training set up for 180s, tone start from 60s and 90s for 10s and shock will be administered during the last 1s of tone.
4. After finish remove mouse carefully and gently from chamber and return it to the home cage.
5. Before we start with another mouse clean the acrylic chamber and floor with 70% of ethanol carefully.

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Day 2. Contextual test

Context testing is performed first, it is usual to do this at the same time of day as the training. Set up computer program and equipment (Acrylic chamber with shock grid) same chamber with conditioning training. Set up light (100 Lux) using (Shodokei).

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1. Mouse handling for 1 minute then placed mouse carefully to the conditioning chamber.
2. Habituation of chamber before actual test for 120s.
3. Start the contextual testing for 180s, without tone and shock.
4. After finish remove mouse carefully and gently from chamber and return it to the home cage.
5. Before we start with another mouse clean the acrylic chamber and floor with 70% of ethanol carefully.

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Day 3. Tone test

Tone testing it is usual to do this at the same time of day as the training and contextual test. Set up computer program and equipment (different chamber with training and also contextual test/ new chamber is use and put wood chip from the mouse home cage). Set up light (100 Lux) using (Shodokei). Set up 65 dB for tone (same dB with training).

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1. Mouse handling for 1 minute then placed mouse carefully to the new chamber.
2. Habituation of chamber before actual test for 120s.
3. Start the tone testing for 180s, with tone from 60s to 120s without shock
4. After finish remove mouse carefully and gently from chamber and return to the home cage.
5. Before we start with another mouse clean the chamber with 70% of ethanol carefully then change the wood chips every time we use different mouse.

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